
Precautions for Unpacking

CX31


◎ In the condition immediately after the CX31 microscope frame is taken out of the package, the microscope comes with various protection measures to prevent deterioration in quality during transport. First remove the protective materials shown in the following illustration.

★ Since the microscope is a precision instrument, handle it with care so as not to apply impact to it.

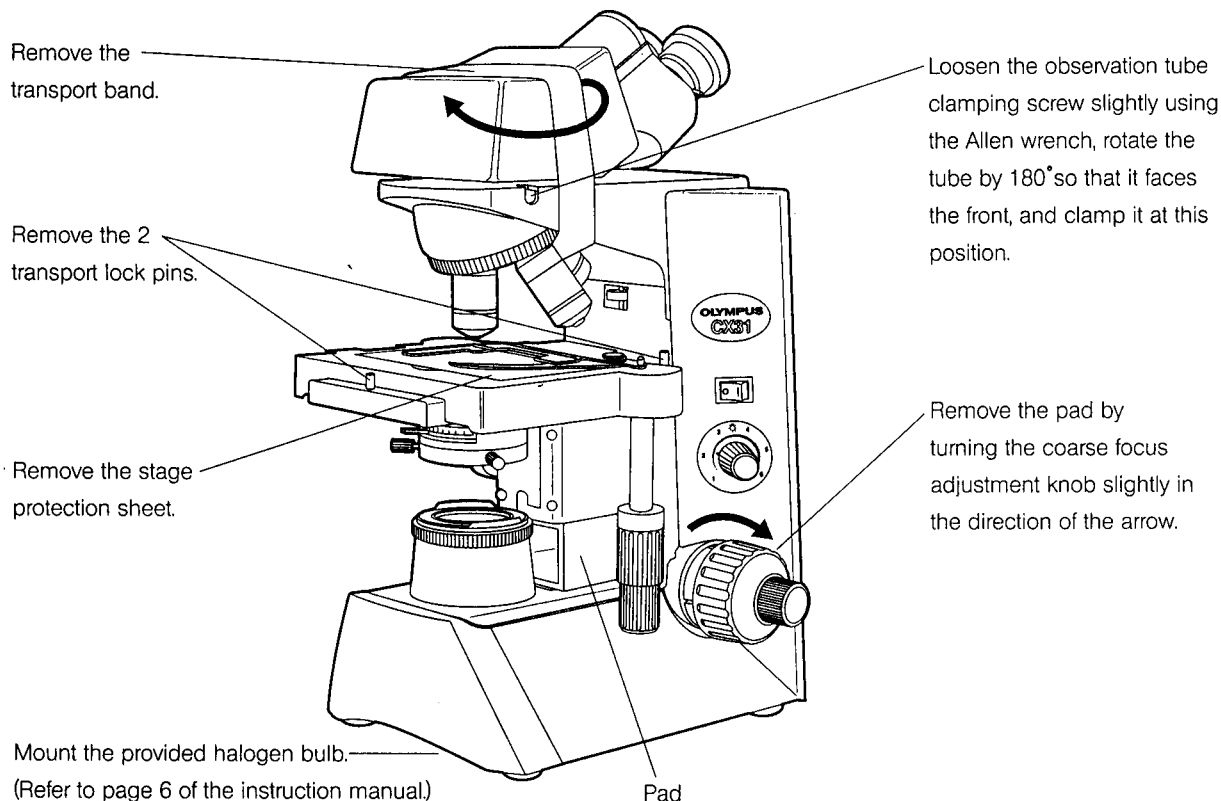◎ The package includes the Instruction Manual, halogen bulb (6V, 30W), Allen wrench and immersion oil (when you purchase a 100X objective) as well as the CX31 microscope frame.

How to Hold the Microscope Correctly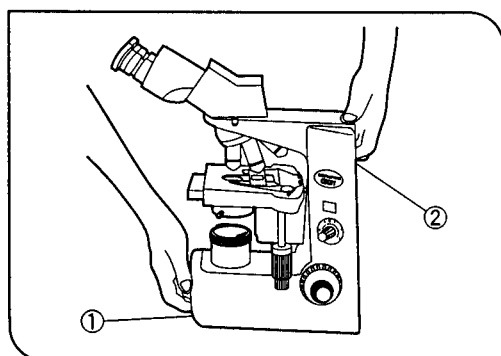

◎ When moving the microscope, carefully carry it with one hand under the base ① and the other hand holding at the recessed handle on the rear of the arm ② as shown in the illustration on the left.

★ Damage to the microscope will occur if you hold it by the stage, X-axis/Y-axis knob, binocular section of the observation tube, etc. Also make sure that specimen, filters, etc. do not fall off.

★ Sliding the microscope on the surface of the table may damage or tear off the rubber feet and/or scratch the table top surface.

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OLYMPUS OPTICAL CO.,LTD.